

Aperio Spectrum User Guide

Please contact us at hudsen@ncl.ac.uk to obtain a user name and password.


After logging in:

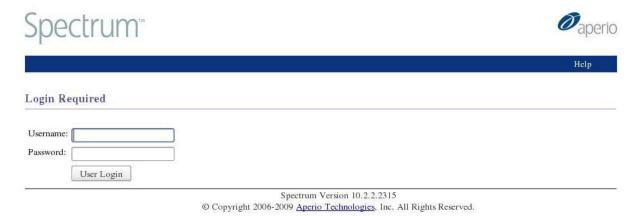

A new window opens. Each set of H&Es is displayed as a folder on the left hand side. By default 20 folders are displayed per page – there may be more than one page of results.

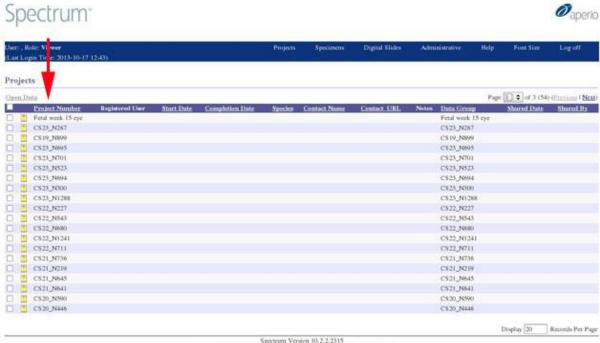

Clicking on the yellow folder icon next to the stage and embryo number brings up thumbnails of all the scanned slides from that embryo.

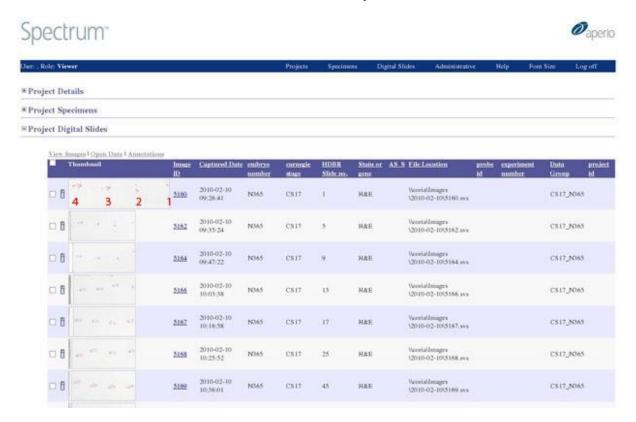



Sections are usually mounted 4 per slide. The first section is furthest away from the frosted end of the slide and appears here on the right hand side of the image (as shown above in red).

Access the scanned slide by clicking on the thumbnail.



By using the tools at the bottom of the page (circled in red above) you can pan across the slide and zoom in up to the full resolution of x20.

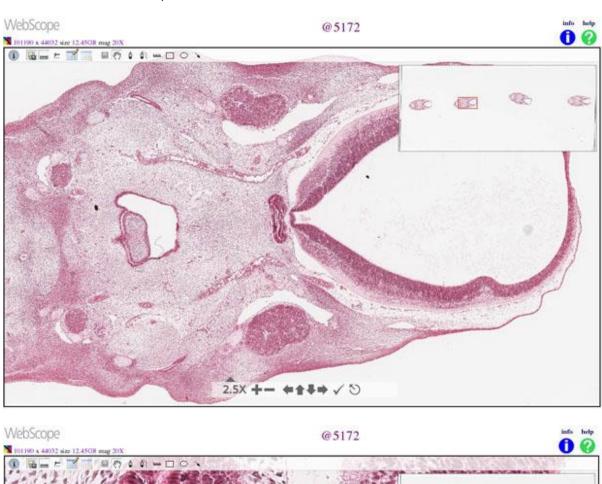

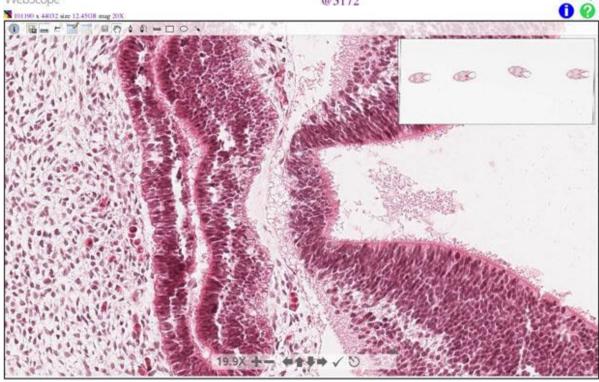

For further information please contact us at husden@ncl.ac.uk.